

Platinum[®] PCR SuperMix High Fidelity

Cat. No. 12532-016

Size: 100 Reactions

Cat. No. 12532-024

Size: 5000 Reactions

**Store at -20°C in a
non-frost-free freezer.**

Description:

Platinum[®] PCR SuperMix High Fidelity provides qualified reagents for high fidelity amplification of DNA templates by polymerase chain reaction (PCR). It is effective over a large range of target sizes—up to 15 kb genomic DNA. The mixture contains anti-*Taq* DNA polymerase antibody, Mg⁺⁺, deoxyribonucleotide triphosphates, recombinant *Taq* DNA polymerase, and *Pyrococcus* species *GB-D* thermostable polymerase. Anti-*Taq* DNA polymerase antibody inhibits polymerase activity, providing an automatic “hot start” (1,2) and permitting room temperature set-up (polymerase activity is restored after a denaturation step in PCR cycling at 94°C). Antibody-mediated hot starts improve PCR specificity and yield (3). *Pyrococcus* species *GB-D* polymerase possesses a proofreading ability by virtue of its 3' to 5' exonuclease activity (3). Mixture of the proofreading enzyme with *Taq* DNA polymerase increases fidelity approximately six times over that of *Taq* DNA polymerase alone.

Platinum[®] PCR SuperMix High Fidelity is supplied at 1.1X concentration to allow approximately 10% of the final reaction volume to be used for the addition of primer and template solutions. Reagents sufficient for 100 or 5000 amplification reactions of 50 µl each are provided.

Component:

100 Rxn Kit

5000 Rxn Kit

Platinum[®] PCR SuperMix High Fidelity*

4 × 1.125 ml

4 × 56.25 ml

*22 U/ml complexed recombinant *Taq* DNA polymerase, *Pyrococcus* species *GB-D* thermostable polymerase, and Platinum[®] *Taq* Antibody; 66 mM Tris-SO₄ (pH 8.9); 19.8 mM (NH₄)₂SO₄; 2.4 mM MgSO₄; 220 µM dNTPs; and stabilizers.

Part. No. 12532

Doc. Rev. 090502

Storage Conditions:

After thawing, Platinum® PCR SuperMix High Fidelity may be stored at 4°C for 3 months or -20°C for 1 year. Storage at 4°C avoids the necessity of thawing the mix before assembling the reaction. There is no detectable decrease in enzyme activity or performance after storage for 3 months at 4°C, or after 15 freeze-thaw cycles.

Recommendations and Guidelines:

- Because PCR is a powerful technique capable of amplifying trace amounts of DNA, take all appropriate precautions to avoid cross-contamination. Ideally, amplification reactions should be assembled in a DNA-free environment.
- Reactions may be assembled either at room temperature or on ice. We have observed no significant difference in reaction efficiency between these setup conditions.
- For multiple reactions, you can prepare a master mix of Platinum® PCR SuperMix High Fidelity and the component(s) common to all reactions.

PCR Protocol:

1. Add the following components in any order to each reaction tube:
 - a. 45 µl Platinum® PCR SuperMix High Fidelity
 - b. Primer solution (200 nM final concentration of each is recommended)*
 - c. Template DNA solution (1–200 ng genomic DNA)*

*A standard 50-µl PCR reaction includes a combined primer and template volume of 5 µl; we have observed no decrease in product yield if the amount of primer and template solution is between 1 µl and 15 µl.
2. Mix contents of tubes and cover with mineral or silicone oil if necessary.
3. Cap tubes and load in thermal cycler.
4. Incubate tubes at 94°C for 30 s to 2 min to completely denature the template and activate the enzyme.
5. Perform 25-35 cycles of PCR amplification as follows:

Denature	94°C for 15-30 s
Anneal	55°C for 15-30 s
Extend	68°C for 1 min per kb

Tips:

- If the PCR efficiency is not optimal, repeat the reaction with different primer concentrations from 100 to 500 nM, in 100 nM increments.
- For longer genomic DNA targets (>15 kb), we recommend adding 1–1.5 U of Platinum® *Taq* DNA Polymerase (Cat. no. 10966-018) to the reaction mix and increasing the extension time as specified (1 min per kb).
- At higher volumes of primer and template, the MgSO_4 concentration in the reaction will drop to suboptimal levels and yield will decrease. For combined primer-template volumes of >15 μl (in solution with 45 μl of Platinum® PCR SuperMix High Fidelity), we recommend adjusting the final MgSO_4 concentration in the reaction to 2.2 mM.

Quality Control:

Platinum® PCR SuperMix High Fidelity is evaluated in a DNA polymerization activity assay that measures the percent of *Taq* DNA polymerase inhibition versus an uninhibited control. A functional assay is also performed. Components are tested for the absence of DNase, RNase and exonuclease activities. Recombinant *Taq* DNA polymerase is tested for the absence of exonuclease, and double- and single-stranded endonuclease activities. The enzyme is >90% homogeneous as determined by SDS-polyacrylamide gel electrophoresis.

References:

1. Chou, Q., Russel, M., Birch, D., Raymond, J., Bloch, W. (1992) *Nucl. Acids Res.*, 20, 1717.
2. Sharkey, D.J., Scalice, E.R., Christy, K.G., Atwood, S.M., Daiss, J.L. (1994) *BioTechnology*, 12, 506.
3. Westfall, B., Sitaraman, K., Solus, J., Hughes, J., Rashtchian, A. (1997) *Focus*®, 19,2, 46.

Limited Use Label License No. 2: PCR Kits

A license under U.S. Patents 4,683,202, 4,683,195 and 4,965,188 or their foreign counterparts, owned by Roche Molecular Systems, Inc. and F. Hoffmann-La Roche, Ltd. ("Roche"), has an up-front fee component and a running-royalty component. The purchase price of this product includes limited, nontransferable rights under the running-royalty component to use only this amount of the product to practice the Polymerase Chain Reaction (PCR) and related processes described in said patents solely for the research and development activities of the purchaser when this product is used in conjunction with a thermal cycler whose use is covered by the up-front fee component. Rights to the up-front fee component must be obtained by the end user in order to have a complete license. These rights under the up-front fee component may be purchased from Applied Biosystems or obtained by purchasing an Authorized Thermal Cycler. No right to perform or offer commercial services of any kind using PCR, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is hereby granted by implication or estoppel. Further information on purchasing licenses to practice the PCR process may be obtained by contacting the Director of Licensing at Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404 or at Roche Molecular Systems, Inc., 1145 Atlantic Avenue, Alameda, California 94501.

Limited License No. 12: SuperMix Products

The purchase of this product conveys to the buyer the non-transferable right to use the product and components of the product in research conducted by the buyer. The buyer cannot sell or otherwise transfer this product or its components to a third party and in particular, no rights are conveyed to the buyer to use the product or its components for commercial purposes. Commercial purposes means any activity for which a party receives consideration and may include, but is not limited to, (1) use of the product or its components in manufacturing, (2) use of the product or its components to provide a service, information or data, (3) use of the product or its components for diagnostic purposes, or (4) resell the product or its components, whether or not such product or its components are resold for use in research. If purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the products with a full refund. For information on purchasing a license to this product for purposes other than research, contact the Director of Licensing, 1600 Faraday Avenue, Carlsbad, California 92008. Phone (760) 603-7200. Fax (760) 602-6500.

Limited Label License No. 14: Platinum[®] products

Licensed to Invitrogen Corporation, under U.S. Patent Nos. 5,338,671; 5,587,287, and foreign equivalents for use in research only.