# **Miniprep - GET Buffer**

### **Solutions/reagents:**

- o overnight culture
- o ice-cold GET buffer

(50 mM glucose (MW 180), 10mM EDTA, 25 mM Tris-HCl pH 8)

- o 0.2M NaOH stored at room temperature
- o ice-cold potassium acetate solution

(3 M potassium acetate, 1.8 M acetic acid, no pH adjustment)

- o 95% / 100% ethanol
- o 70% EtOH
- o TE buffer
- o distilled water
- o PCA solution

((optional)50 parts phenol, 49 parts chloroform, and 1 part amyl-alcohol)

- o SDS
- lysozyme (optional)

### **Steps:**

- 1. Measure out 2 ml of overnight culture into a sterile 2-ml microcentrifuge tube.
- 2. Centrifuge at maximum speed for 1 min at room temperature, gently aspirate out the supernatant and discard it.
- 3. Add 100 µl of ice-cold GET buffer.

Resuspend the pellet by vortexing/by shaking vigorously.

### (Optional)

Add 10 mg of lysozyme.

Store at room temperature for 30 mins.

This step is essential for lysing gram-positive cells.

4. Measure out  $200~\mu l$  of 0.2M NaOH into a sterile 2-ml microcentrifuge tube.

Add 2 mg of SDS.

Vortex the mixture for a few secs.

5. Add alkaline SDS solution to cell suspension.

Close the tube tightly and gently mix the contents by inverting the tube.

DO NOT VORTEX! The solution should become clear.

6. Add 150 μl of ice-cold potassium acetate solution.

Close the tube tightly and gently mix the contents by inverting the tube.

DO NOT VORTEX! A precipitate should form.

- 7. Store the tube on ice for 3 5 mins.
- 8. Centrifuge at maximum speed for 10 mins at room temperature and aspirate out 400 µl of top layer.

Transfer top aqueous layer into a sterile 2-ml microcentrifuge tube.

Discard bottom layer.

DO NOT PICK UP ANY PRECIPITATE!!!

#### (Optional)

Add 400 µl of PCA solution.

Close the tube tightly and gently mix the contents by inverting the tube.

Centrifuge at maximum speed for 3 mins at room temperature and aspirate out the top layer.

Transfer top aqueous layer into a sterile 2-ml microcentrifuge tube.

Discard bottom layer.

This helps remove any residual proteins.

9. Add **900 μl** of 95% / 100% ethanol.

This is to precipitate the plasmid DNA.

10. Store at -80°C for 30 mins.

Use the -80°C freezer.

11. Centrifuge at maximum speed for 10 mins at room temperature, gently aspirate out the supernatant and discard it.

## **Equipment:**

- o Centrifuge
- o Sterile 2-ml microcentrifuge tubes

- Store at **room temperature** for **3 mins**.

  13. Centrifuge at maximum speed for **3 mins** at **room temperature**, gently aspirate out the supernatant and discard it.
- 13. Centrifuge at maximum speed for 3 mins at room temperature, gently aspirate out the supernatant and discard in Make sure the pellet is toward the outside.
- 14. Dry the pellet in air for 10 15 mins.

12. Add 1 ml of 70% EtOH.

15. Make sure the pellet is completely dry before this step.

Option 1: Add  $20 \mu l$  of TE buffer.

(or

Option 2: Add 20 µl of distilled water.

Resuspend the pellet by vortexing/by shaking vigorously.

The DNA will contain RNA contamination, which can be removed by resuspending in TE with RNAse.