

## Is This Data Right?

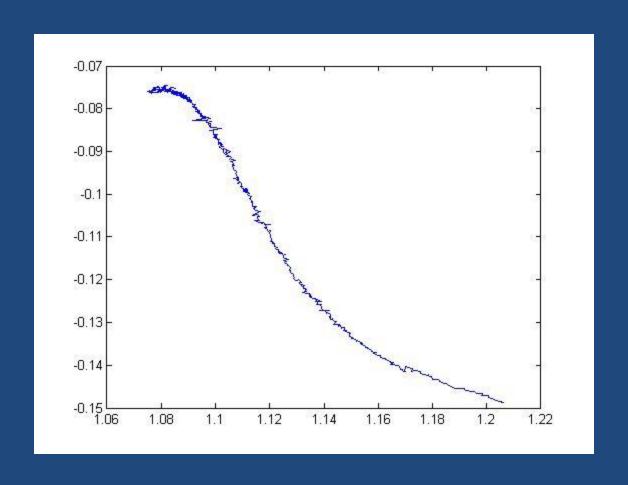

## **Processed Data**

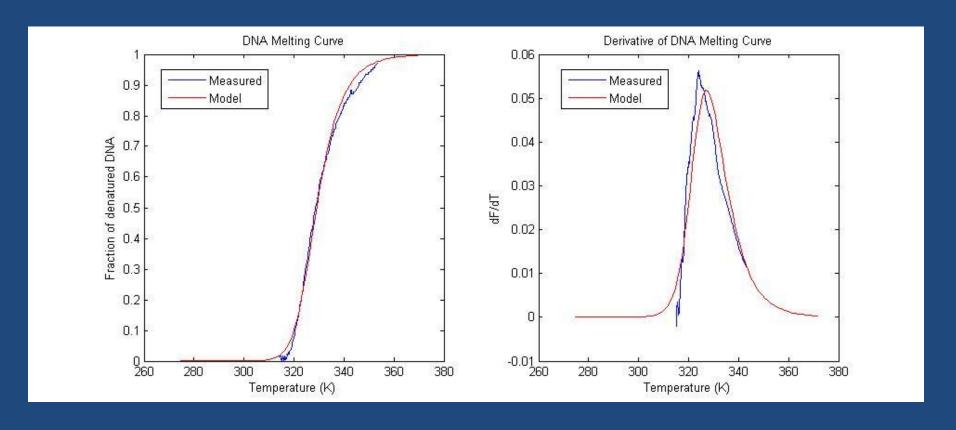

## Analysis

- Filter
  - Median filter to remove impulse noise
    - Replace each sample with the median of it and its neighbors
    - Matlab hint: help medfilt1
  - FIR low-pass filter (averaging) to remove random noise
    - 60 Hz comes through vigorously in the derivative
    - Working on 60 Hz filter in VI
    - Matlab hint: help conv
      - Be careful at the edges conv sets initial conditions to zero
    - Filter design: help fir1
- Transform
  - Range of V<sub>f</sub>
    - Assumption?
  - Convert V<sub>RTD</sub> to temperature
- Take finite difference (derivative)
  - $-\Delta F/\Delta T$
  - Is your data a function?
  - Matlab hint: help diff
- Fit to model
  - Matlab hint: help lsqcurevefit
- Compute melting temperature
  - Different methods?
    - *f=0.5*
    - Maximum df/dt
    - From fit

## Lab Stuff

- Default salt concentration of samples is 100mM
- Dispose of DNA properly
- Please do not drink the DNA
  - Limited supply going astonishingly fast
- The glass cuvettes are expensive be careful
- Clean cuvette with DI water
  - It's okay if the glass cuvette has a few droplets of water in it when you change samples