Reagents

diH₂O

50% 100 mM Ammonium Bicarbonate (AB) / 50% Acetonitrile (ACN)

100% 100 mM AB

100% ACN

10mM Dithiothreitol (DTT) in 100 mM AB

55mM Iodoacetamide (IAA) in 100 mM AB

100% 25 mM AB

Trypsin Digestion Solution (make fresh): Thaw a 50 μL vial of 0.1 ug/ μL trypsin, and

add 200 μL of 100 mM AB, 4 μl 0.5 M CaCl₂,

and 146 $\mu L \, di H_2 O$

Digestion Buffer: 5.0 mL of 100 mM AB, 100 µL of 0.5 M CaCl₂, 4.90 mL H₂O

Pre-Digestion

Solutions removed after each incubation unless otherwise noted

(2x) Add diH₂O wash of whole gel for 15 min

Cut out gel bands and put into vials

(2x) Add 100ul of 50% 100 mM AB / 50% ACN and incubate for 30 min on shaker

Put sample in Speed-Vac to remove excess solution

Add 100 µL of 10 mM DTT and incubate for 45 min @ 56°C

Add 100 µL of 55 mM IAA cover with foil and incubate for 30 min on shaker

Add 500 µL of 50% 100 mM AB / 50% ACN and incubate for 15 min on shaker

Put sample in Speed-Vac to remove excess solution

Digestion

Add 20 μ L of Trypsin Digestion solution and incubate for 15 min @ 4°C If gel plug isn't completely re-swelled add an additional 10 μ L of Trypsin solution and repeat the incubation until the gel plug won't absorb any more solution then remove any excess solution

Add 60 µL of Digestion Buffer and incubate @ 37°C overnight, more digestion buffer can be added to cover the gel slices, don't remove solution after incubation

Short Extraction Method

Add 3 µL of neat FA added to the above solution, bath-sonicate for 15min at 37°C, pull off and save solution for mass spec analysis

Long Extraction Method

Pull off solution from the overnight digestion into a separate vial

Add 50 μ L of 25 mM AB gel pieces, bath-sonicate for 15 min @ 37°C, don't remove solution after incubation

Add 50 µL of 50% ACN to the above solution, bath-sonicate for 15 min @ 37°C, remove supernatant after incubation and combine it with the solution from the overnight digestion

Concentrate the extracts to approx. 10 μL in a Speed-Vac (don't let them go dry) and save for mass spec analysis